

# BIOCLEAR™

## ANTI-FOG SPRAY

For Autoclavable Goggles

# BAFS

### About

To address any fogging issues with our autoclavable goggles an application of BioClean Anti-fog Goggle Spray is recommended. The procedure for the successful application of the spray is as follows:

1. Autoclave the BioClean autoclavable goggles as usual
2. Just before donning, apply one or two sprays to the INSIDE of the lens
3. Using a sterile cleanroom wipe, dry the excess liquid from the inside surface of the lens with a spreading motion, covering the whole of the inside of the lens
4. Ensure the treated side of the lens is dry after wiping
5. Repeat the process after every autoclave cycle

### Specifications

**MATERIAL:**  
Non-hazardous ingredients

**LENGTH:**  
60ml

**COLOUR:**  
Colourless

**STERILIZATION:**  
Gamma irradiation

### Features

- Eliminates fogging
- Airless pump action spray
- Alcohol free
- Quick and easy application
- Suitable for all BioClean autoclavable goggles

**ELIMINATES FOGGING**

**AIRLESS PUMP ACTION SPRAY**

**ALCOHOL FREE**

**QUICK AND EASY APPLICATION**





## Quality Standards

- Manufactured in a facility holding ISO 9001:2008 certifications

## Sterilization

Method: Gamma irradiation  
 Minimum Dose: 25kGy  
 Sterility Assurance Level: 10<sup>-6</sup>

## Shelf Life & Storage

Three (3) years from date of manufacture. Store in a dry, cool place (<40°C) away from direct sunlight and fluorescent light.

## TO ORDER

Re-order code	Size
BAFS-601	60ml

### Packing

One bottle per sealed inner PE bag; 10 inner bags per sealed outer PE bag; 10 outer bags per lined carton (100 bottles)

## CONTACT

### Europe

#### (Headquarters)

Nitritex Ltd.  
 Minton Enterprise Park,  
 Oaks Drive,  
 Newmarket, Suffolk,  
 CB8 7YY,  
 United Kingdom  
 T: +44 (0) 1638 663338  
 F: +44 (0) 1638 668890  
 E: [europa@bioclean.com](mailto:europa@bioclean.com)

### Other Regions

Asia:  
[malaysia@bioclean.com](mailto:malaysia@bioclean.com)  
 Americas:  
[americas@bioclean.com](mailto:americas@bioclean.com)



BioClean™ is a brand of Nitritex Ltd.  
 For more information, please visit [www.bioclean.com](http://www.bioclean.com)

BAFS/1604/PDS3