

BIOCLEAN™

PHARMA COVERS

Equipment & Glassware Covers

BPC

About

BioClean Pharma Covers™ provide protection from environmental impurities during storage, use and transportation. They also minimize the need for autoclave tape during wrapping and save time when prepping glassware, tubing etc for autoclaving. Easy removal minimizes particulate generation and facilitates aseptic technique. Available in diameters 5cm to 229cm, bespoke sizing available on request.

Specifications

MATERIAL:

Tyvek® material and high quality elastic

COLOUR:

White

Features


- Breathable microbial barrier
- Rubber elastic provides secure seal
- Lint-free, moisture and puncture resistant
- Saves time when prepping
- Form fitted
- Fast drying time
- Easy to apply and remove
- Autoclavable at 121°C to 127°C

LINT-FREE

MOISTURE AND PUNCTURE RESISTANT

PROVIDES SECURE SEAL

AUTOCLAVABLE

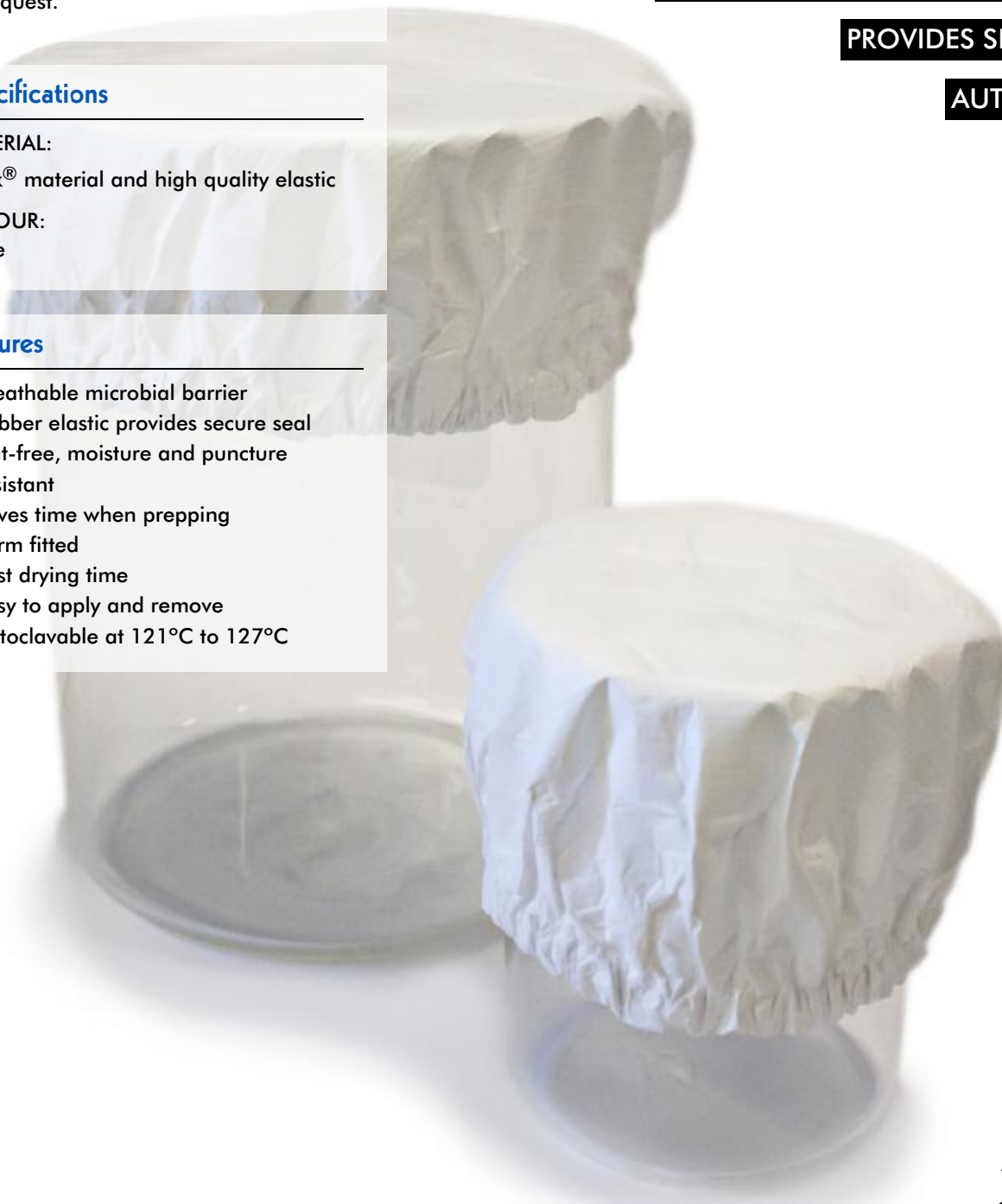

Quality Standards

- Manufactured in a facility holding ISO 9001: 2008 certification

Shelf Life & Storage

Three (3) years from date of manufacture. Store in a dry, cool place (<40°C) away from direct sunlight and fluorescent light.

Test Results

Typical Helmke Drum Test Results			
size(μm)	Average counts per minute		
	BPC-008	BPC-041	BPC-061
≥0.3	104	184	223
≥0.5	93	155	208
≥5.0	36	49	66

Endotoxin Test Results	
Method, USP <85> (Kinetic Turbidimetric Method)	<20 EU/Sample
Bacterial Filtration Efficiency (BFE)	
Method, ASTM F2101	98.8%



TO ORDER

Re-order code	Size
BPC-005	5cm (2.0") ∅
BPC-008	8cm (3.1") ∅
BPC-010	10cm (3.9") ∅
BPC-013	13cm (5.1") ∅
BPC-018	18cm (7.0") ∅
BPC-023	23cm (9.0") ∅
BPC-030	30cm (11.8") ∅
BPC-041	41cm (16.1") ∅
BPC-051	51cm (20.0") ∅
BPC-061	61cm (24.0") ∅
BPC-076	76cm (29.9") ∅
BPC-091	91cm (35.8") ∅
BPC-122	122cm (48.0") ∅
BPC-152	152cm (59.8") ∅
BPC-183	183cm (72.0") ∅
BPC-229	229cm (90.0") ∅

Packing

BPC-005/BPC-008/BPC-010/BPC-013/BPC-018/BPC-023
/BPC-030/BPC-041/BPC-051/BPC-061:

50 pieces per sealed inner PE bag; 10 inner bags per sealed outer PE bag; One outer bag per lined carton (500 pieces)

BPC-076/BPC-091/BPC-122:

10 pieces per sealed inner PE bag; 10 inner bags per sealed outer PE bag; One outer bag per lined carton (100 pieces)

BPC-152/BPC-183/BPC-229:

Two pieces per sealed inner PE bag; 10 inner bags per sealed outer PE bag; One outer bag per lined carton (20 pieces)

CONTACT

Europe

(Headquarters)

Nitritex Ltd.
Minton Enterprise Park,
Oaks Drive,
Newmarket, Suffolk,
CB8 7YY,
United Kingdom
T: +44 (0) 1638 663338
F: +44 (0) 1638 668890
E: europa@bioclean.com

Other Regions

Asia:

malaysia@bioclean.com

Americas:

americas@bioclean.com



BioClean™ is a brand of Nitritex Ltd.

For more information, please visit www.bioclean.com

BPC/1611/PDS26